

If you want to use the Metal frame with membrane I suggest to you to use the Plexiglas FrameSupport (code: 11532325)



This support (transparent with Leica logo) has been included in the delivery system.

**Plexiglass FrameSupport contains a lifted surface (clear area) that fits exactly into the cavity or the front side of the FrameSlide.**

It must be placed exactly at the raised part of the "cavity" of the frame with membrane that contains the Leica logo printed. The section at this point is collected on the flat surface of the frame (the one without Leica logo)

Plexiglas FrameSupport with this momentary adhesion of the metal frame reduces the cooling rate by facilitating the adhesion of frozen sections.

After adhesion of the section to the membrane of metal frame we can remove the metal frame with the membrane from the Plexiglas FrameSupport and we can proceed with the standards staining